

NamaZol RNA Extraction Body Fluids

1

**NamaZol (750µl)
Sample (250µl)**

**Incubate:
RT
5-15 min**

2

**Chloroform (200µl)
Shake for 15 sec**

**Incubate:
RT/2-5 min
Cen:
15 min/12000 rpm/ 4 °C**

3

**Transfer upper phase to new tube
Isopropanol (500 µl)
Nucleic acid carrier (1-10 µl)**

**Incubate:
RT/5-10 min
Cen:
10 min/12000 rpm/ 4 °C**

4

Ethanol 75% (1ml)

**Cent:
5min
8000 rpm/4 °C**

5

**Remove Ethanol & air-drying
DEPC-trated water (20-100µl)**

**Incubate:
55 °C
10-15 min**